

ROTARY EVAPORATOR



STRIKE[®]100

STRIKE100 is a large-capacity rotary evaporator designed for industrial plants. It meets all the process needs by using Borosilicate glass and PTFE permitting aggressive substances treatment.

The unit works under vacuum or atmosphere conditions. The rotation of the 50 or 100 lt evaporation flask in the thermostatic bath results in a forced convection and homogenous distribution of the product, thus preventing sedimentation. Moreover, it ensures high evaporating exchange surface. The sealing system guarantees the perfect vacuum tightness and maintenance-free operation. Possibility to customize glassware as for Clients needs.

APPLICATIONS

- Distillation (continuous and discontinuous)
- Solvent recycling
- Reflux process reactions
- Component drying
- Fine chemical synthesis
- Crystallization

FIELDS

- Pharmaceutical
- Chemical
- Cosmetic
- Herb extraction
- Petrochemical



TOTAL SAFETY SYSTEM

The industrial safety rotary evaporators are able to safeguard the user from any risk during work. The thermostatic bath is covered by a transparent shield providing the operator with utmost protection from glass breaking and spray at high temperature. The thermostatic bath and the protective shield are lifted by means of an hydraulic system electronically controlled. An automatic system lowers the thermostatic bath when a power failure, door opening and overheating occur. It is also inserted on glassware a brake disc in order to avoid sudden overpressure.

VACUUM SEALING SYSTEM

A great feature of these rotary evaporators is its sealing system made of self-lubricating material which allows a perfect vacuum-tight. This sealing system does not require any maintenance.



HEATING BATH

The heating bath is lifted by means of a hydraulic system which, in case of power failure or doors opening, goes automatically down to avoid fluid overheating and the process will immediately stop. Electric heater enhanced to 8Kw to ensure a continuous evaporation of the product even in the worst conditions.

DISPLAY

The process operating conditions can be set and displayed: any value error or anomalous temperature will block the electric power. For STRIKE100 ATEX version is a "LCD display anti-reflectio coating glass panel!". For version STRIKE100 is a graphical color display 7" touch-screen

ROTARY EVAPORATOR



STRIKE[®] 100

ELECTRONIC CONTROLLER


Strike100 e STRIKE100ATEX are fully controlled by a PLC (Programmable Logic Controller) which allows for controlling the parameters as well as setting them directly by means of the keys.

Client could acquire on an external PC through serial communication all the process parameters through special software for data acquisition and management. Client could remotely control the instrument from an external PC via serial communication. Client could save and recall up to 20 methods or "recipes" work that the customer can customize over time (NOT ON ATEX VERSION). In addition, the instrument has control algorithms that allow you to finish the process of evaporation / distillation or separation of solvents automatically by reading the temperature of the vapors.



PARAMETERS MEASUREMENT, VISUAL DISPLAY AND RECORDING OF THE FOLLOWING:

- Bath pre-set and real temperature
- Vapor pre-set and real temperature
- Pre-set and real rotation speed
- Pre-set and real vacuum reading and setting

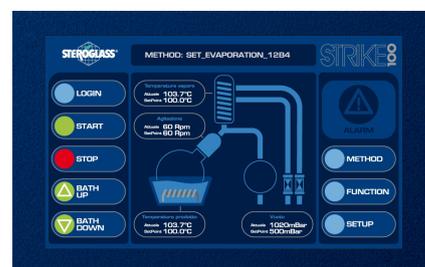

HOW TO ORDER

Code	STRIKE 100	Code	STRIKE 100 ATEX *
	100 lt version		100 lt version
SQFU069840	100lt reflux glassware 1,43m ² 2x20lt	SQFU069848	100lt reflux glassware 1,43m ² 2x20lt
SQFU069841	100lt reflux glassware 1,43m ² 1x20lt	SQFU069849	100lt reflux glassware 1,43m ² 1x20lt
SQFU069842	100lt descending glassware 100lt 1,43m ² 2x20lt	SQFU069850	100lt descending glassware 100lt 1,43m ² 2x20lt
SQFU069843	100lt descending glassware 100lt 1,43m ² 1x20lt	SQFU069851	100lt descending glassware 1,43m ² 1x20lt
	50 lt version		50 lt version
SQFU069844	50lt reflux glassware 0,6m ² 2x20lt	SQFU069852	50lt reflux glassware 0,6m ² 2x20lt
SQFU069845	50lt reflux glassware 0,6m ² 1x20lt	SQFU069853	50lt reflux glassware 0,6m ² 1x20lt
SQFU069846	50lt descending glassware 0,6m ² 2x20lt	SQFU069854	50lt descending glassware 0,6m ² 2x20lt
SQFU069847	50lt descending glassware 0,6m ² 1x20lt	SQFU069855	50lt descending glassware 0,6m ² 1x20lt

*Strike 100 Atex version is in compliance with the Atex European Directives that regulate explosion-proof manufacturing materials. Strike 100 Atex is T3 or T4 certificated. T3 or T4 protectionlevel can be chose according to the applications needed.

ROTARY EVAPORATOR



STRIKE®

Technical features

Thermic Bath		Interface	
Working Range °C	[0; 150]	PLC controlled	YES
Accuracy °C	+/- 2	Export data interface	YES
Resolution °C	0.1	Remote control interface	YES
Set point	YES (1°C step)	Alarms	
Type of control	PID	Thermal bath overheat	YES
Maximum oscillation during checkout °C	+/- 3	Bath temperature exceeding set point	YES
Power resistor KW	8	Soft start rotation	YES
Heater power density	3W/cm2	Rotation speed monitoring	YES
Over-temperature protection	YES	Security status (rotation, lifting, heating=off)	YES
Temperature Sensor	Three-wire PT100	Evaporation flask protection opening	YES
Level sensor	YES	General	
Filling Level	Yes -Manual	100Lt evaporation flask	YES
Vapour Temperature		50Lt evaporation flask	YES
Working Range °C	[0; 150]	20L evaporation flask	YES
Accuracy °C	+/- 2	20Lt collecting flask	2
Resolution °C	0.1	cooler exchange surface (mq)	0.6/1.4/custom
Temperature sensor	Three-wire PT100	Over-pressure Protection	YES
Rotation		Protection with rupture disk	YES
Working Range Rpm	[5; 100]	Quick closing	YES
Accuracy Rpm	+/- 2	ATEX certification	YES
Rotation speed reading	YES	Dimensions (WxDxH mm)	1830x890x2110
Resolution Rpm	1	Operation	
Engine 0,18KW, 380Vac	YES	vapours cycle maximum temperature	YES
Vacuum Control		vapours cycle minimum temperature	YES
Working Range mbar	[0; 1000]	vapours cycle Temp range	YES
Accuracy mbar	±5mbar	receipt/methods in memory	YES
Resolution mbar	1	recipes customization	YES
Absolute vacuum reading	YES	Data acquisition on PC	YES
Display	Digital	Remote Control	YES
Vacuum regulation	YES	lowering bath in case of alarm	YES
Lifting		Optional	
Lifting control ON/OFF	YES	Plastic coated glassware	
Hydraulic lifting	YES		



STEROGLASS S.r.l.

Strada Romano di Sopra 2/C, 06132 San Martino in Campo (PG), Italia
Tel. +3907560909-1 fax. +390756090950 e-mail. info@steroglass.it

ROTARY EVAPORATOR



STRIKE®

STEROGLASS®

COMPANY WITH
QUALITY SYSTEM
CERTIFIED BY DNV
= ISO 9001 =

STEROGLASS s.r.l.
Via Romano di Sopra, 2/C

06132 - S.Martino in Campo
PERUGIA - ITALY
Tel. +39 075 60 90 91
Fax +39 075 60 90 950

http: www.steroglass.it
e-mail: info@steroglass.it

Authorized Dealer

